

Glycoprotein staining- PAS Schiff's reagent

- 1- 200 ml 40% Methanol and 7% acetic acid – 1hr RT
- 2- (800 ml) , used 100 ml 7.5% acetic acid- 1 hr RT
- 3- 200 ml 1% periodic acid 1 hr (4 C, dark)
- 4- Wash in 7.5% Acetic acid for 10 min 6X
- 5- Add 50-100 ml Schiff's reagent- 1 hr (4 C in the dark)
- 6- (400 ml) Wash in 100 ml 0.5% sodium metabisulphate
- 7- Rinse in d-water